IMMB - 0 9 - 02- 0 5 -22- EN -IT

Startup Guide

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Precision Balances

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-
- PM C32

To view a complete User Manual go to the website or scan QR code:

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
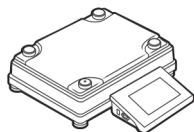
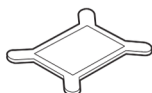
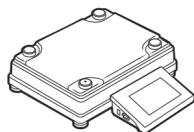
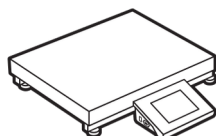
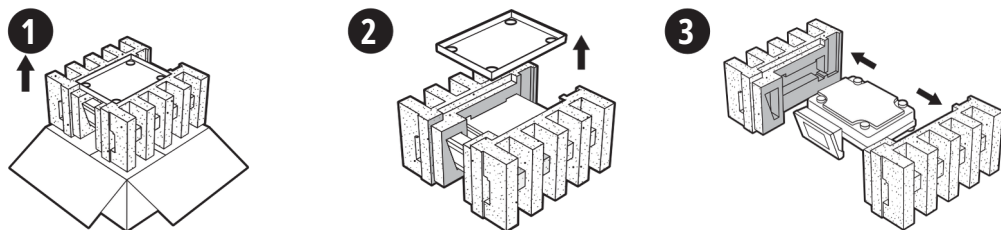
PM C32 (max: 10-50 kg)

PM C32 (max ≥ 60 kg)

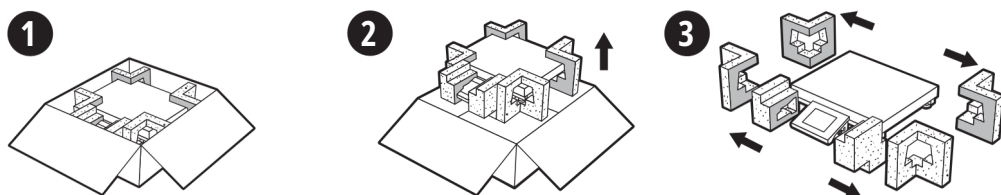
Keep the packaging should a warranty claim or service be required.

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3. RECOMMENDATIONS AND WARNINGS

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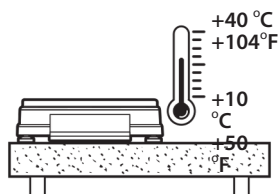Operate the device in a room where the temperature ranges between

10–40 °C (50–104 °F) and where the relative humidity is below 80%.

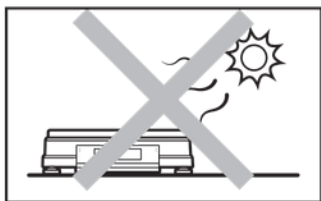
Temperature must change slowly and gradually.

Using the anti-vibration table is recommended in order to enable the most stable and repeatable measurement.

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

Place the balance away from heat sources. Avoid exposing the balance to the sunlight.

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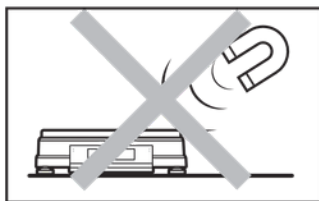
Avoid exposing the balance to a magnetic field. Do not weigh magnetic substances.

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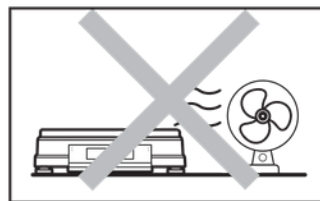
Avoid air drafts and air movements at the workstation.

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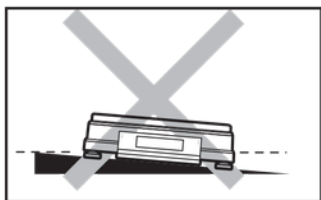
Make sure that the balance is placed on an even surface.

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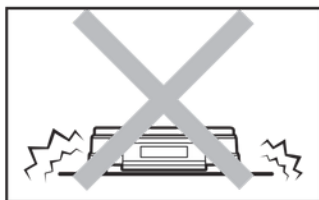Do not place the balance on an unstable surface exposed to shocks and vibrations.

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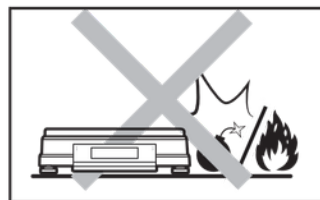Not intended for usage within EX zones. Not intended for weighing explosive or flammable substances.

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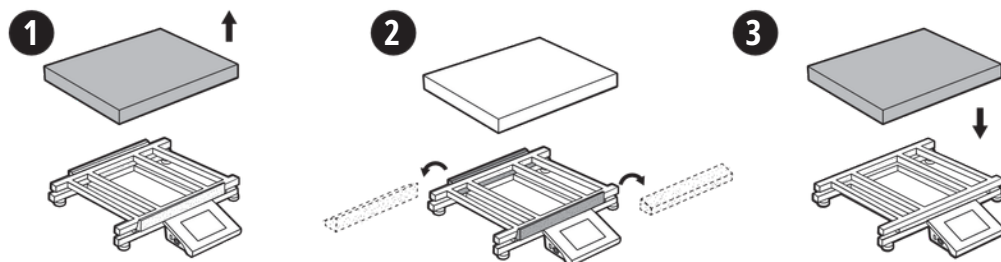
4. ACTIVITIES TO BE DONE PRIOR TO OPERATION

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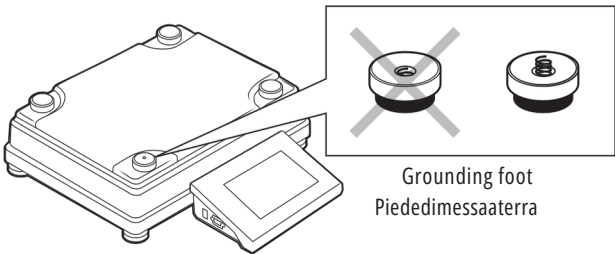
PM C32 (max: 10-50 kg)


Check whether the ground-
ing spring takes its intended
location. Make sure that the
grounding spring juts slightly
outofthe hole.

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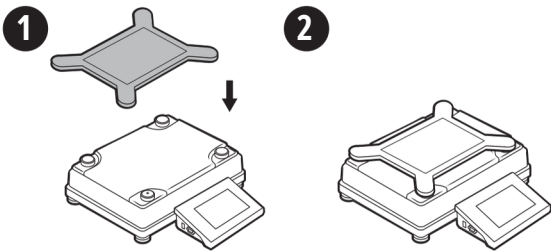
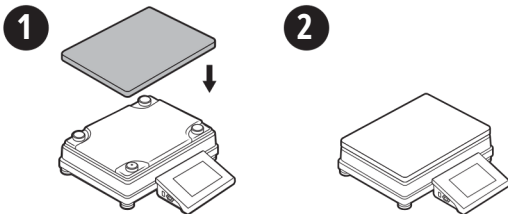
PM C32 (max ≥ 60 kg)

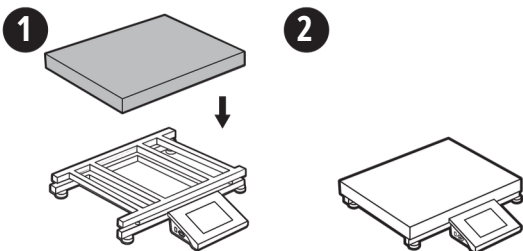6. PREPARING FOR WORK

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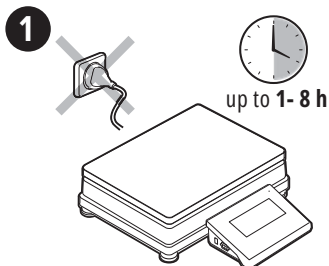
Place the device at the workstation, wait until it reaches room temperature.

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

Prior to switching the device on, let it reach room temperature (estimated stabilisation time: 1-8 hours).

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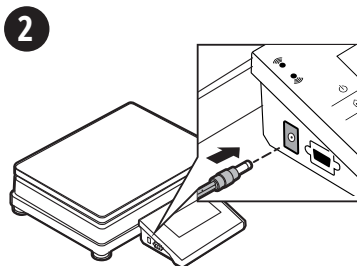Connect the power adapter to DC connector.

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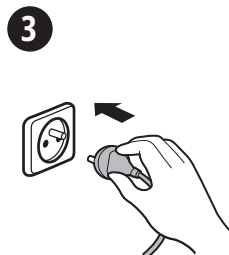
Connect the power adapter to the mains.

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Livellamento

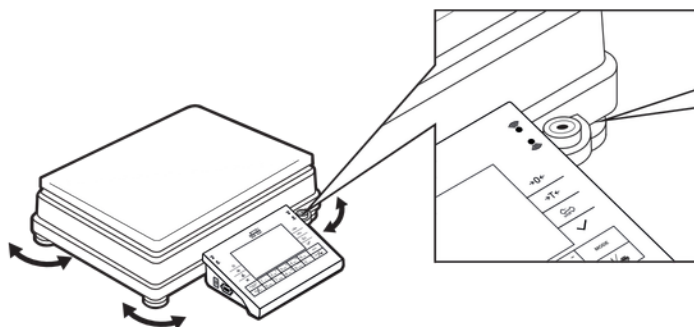
PM C32 (max: 10 - 50 kg)

To level the balance, turn its feet and keep turning the feet until an air bubble takes the central position.

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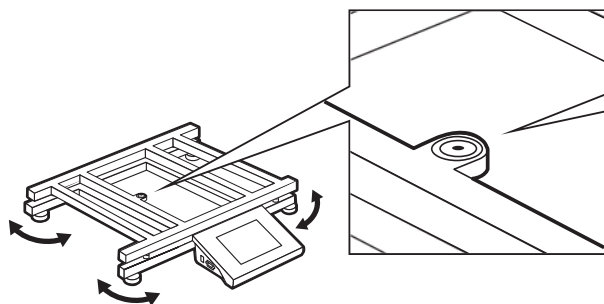
6.3. Switching the balance on

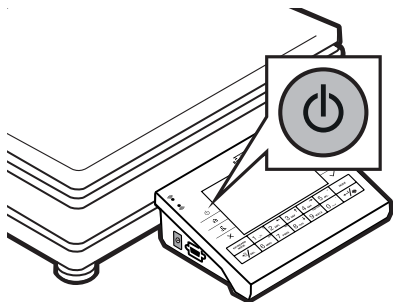Accendere la bilancia

To switch the device on, press ON / OFF key.

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

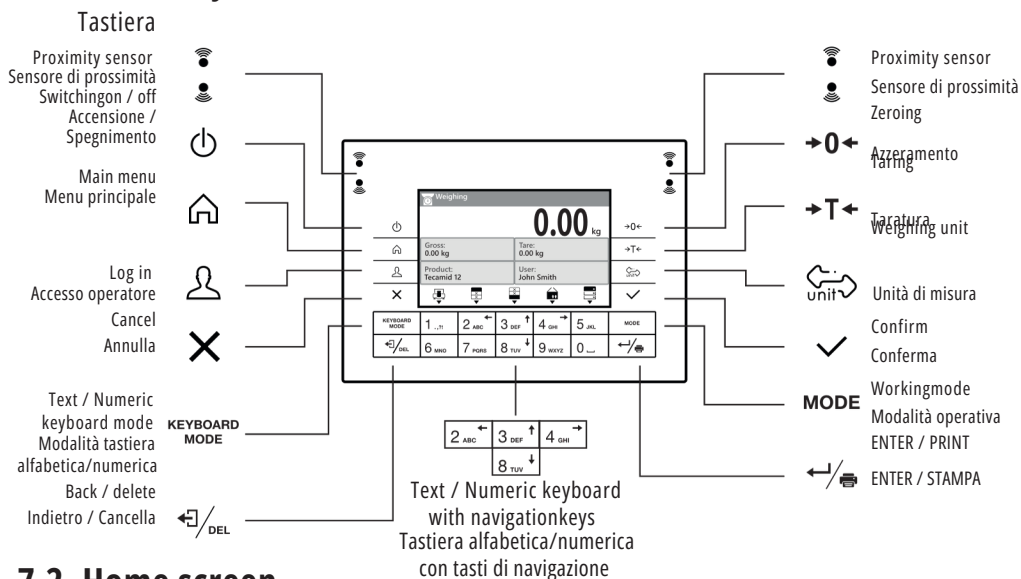
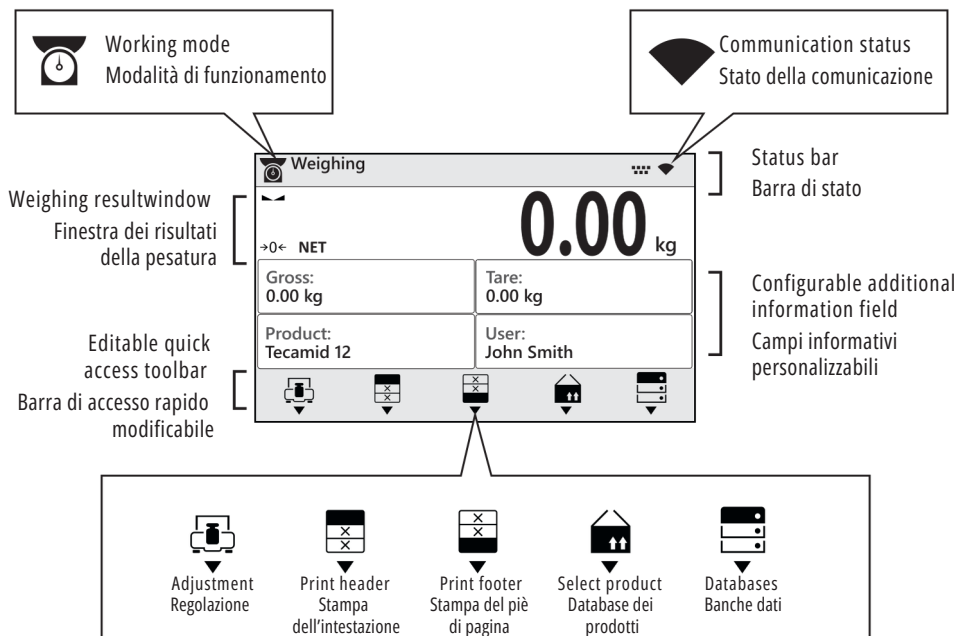
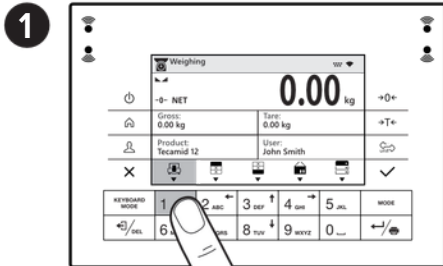
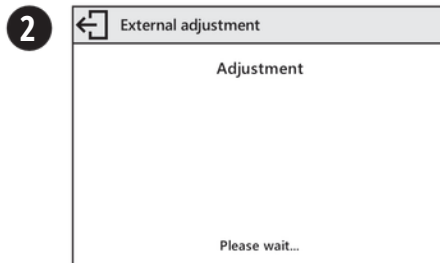
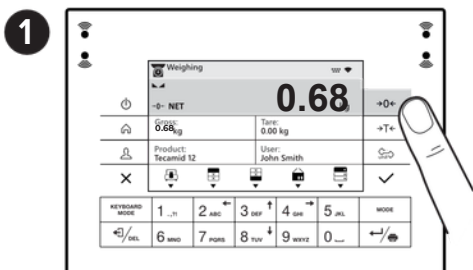
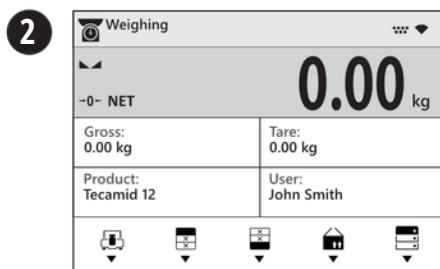
8.3. Taring

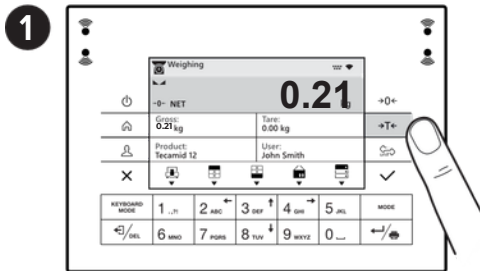Taratura

With loaded weighing pan: upon stabilization of the weighing result, press **+T** key.

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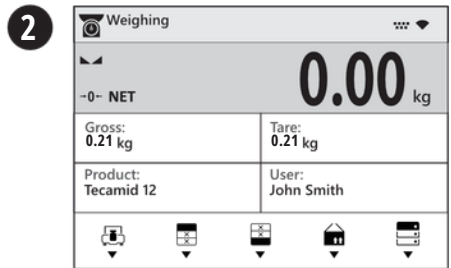
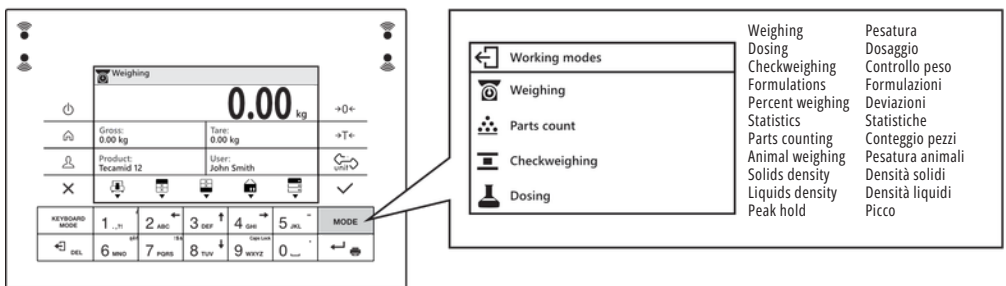
8.5. Units selection

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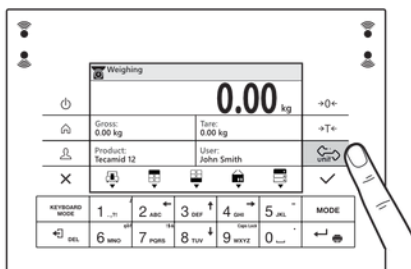Press the **UNIT** key to change the weight unit.

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The balance enables selecting the following weighing units:

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9. SETTINGS

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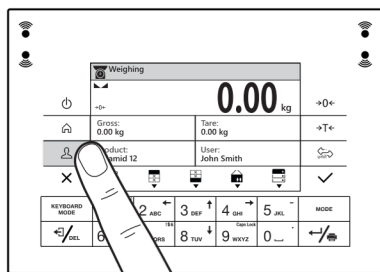
**Some balance settings are accessible for Administrator exclusively.
Prior balance parameters setup, login as the Administrator.**

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
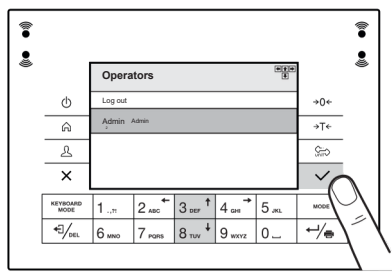
2

Users list is displayed. Using  keys,
select "Admin" option and press  key
to confirm.

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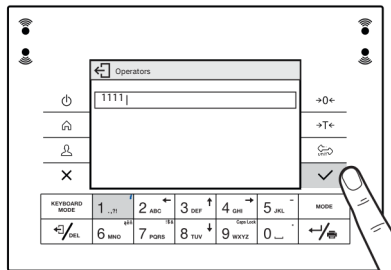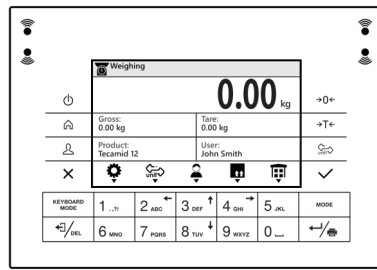
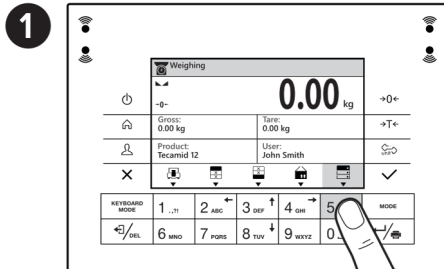
9.2. User settings

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
Prior balance users setup, log in as Administrator (point 9.1.).

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
Press the key under the  icon to enter databases.

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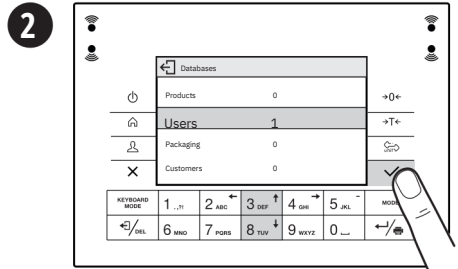
Using  keys, go to "Users" and press  key to confirm.

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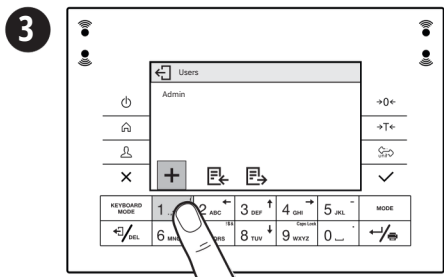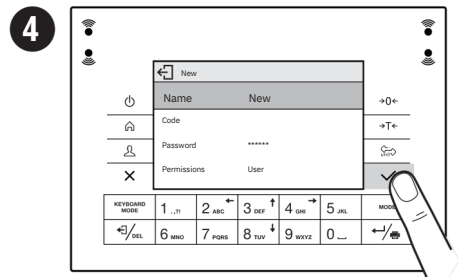
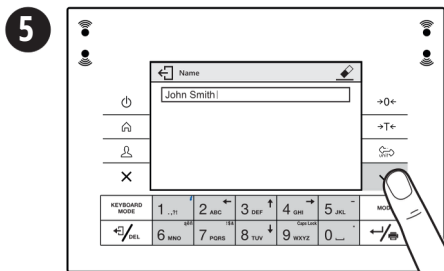
Using keyboard enter user name and press  key to confirm.

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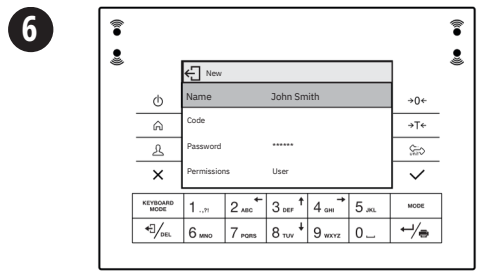7

Following the above procedure, set remaining parameters:

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Set respective access level (user, advanced, admin)

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New	
Name	John Smith
Code	
Password	*****
Permissions	User

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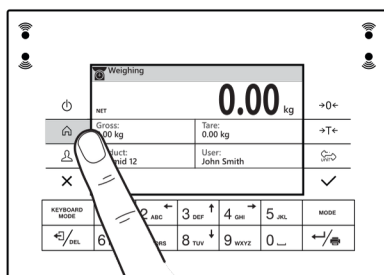
Prior balance proximity sensors setup, login as Administrator (point 9.1.).

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You can assign the proximity sensors with different functions for each working mode (the sensors will trigger different operation for each working mode). The diagram presents sensor setup for the "Weighing" mode.

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

Premere il pulsante .

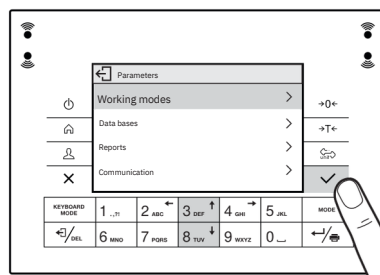2


Using   keys, go to "Working Modes" and press  key to confirm.

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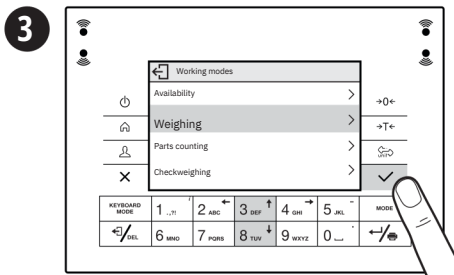
Using keys select "Weighing" mode, and press key to confirm.

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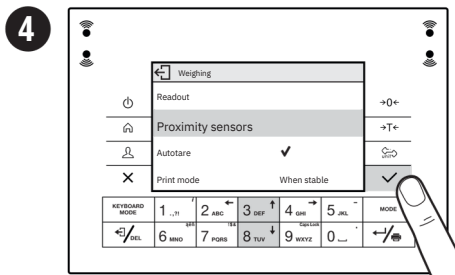
Using keys select "Proximity sensors", and press key to confirm.

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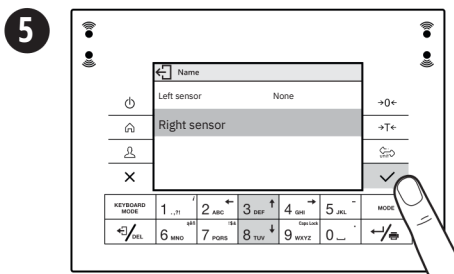Using keys select the appropriate sensor (right or left), and press key to confirm.

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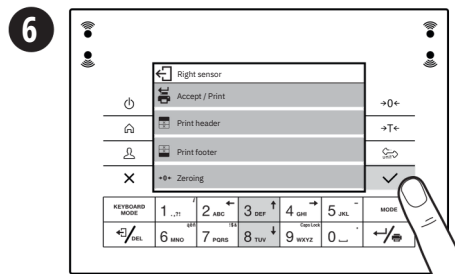Using keys set function that is to be assigned to a selected sensor, and press key to confirm.

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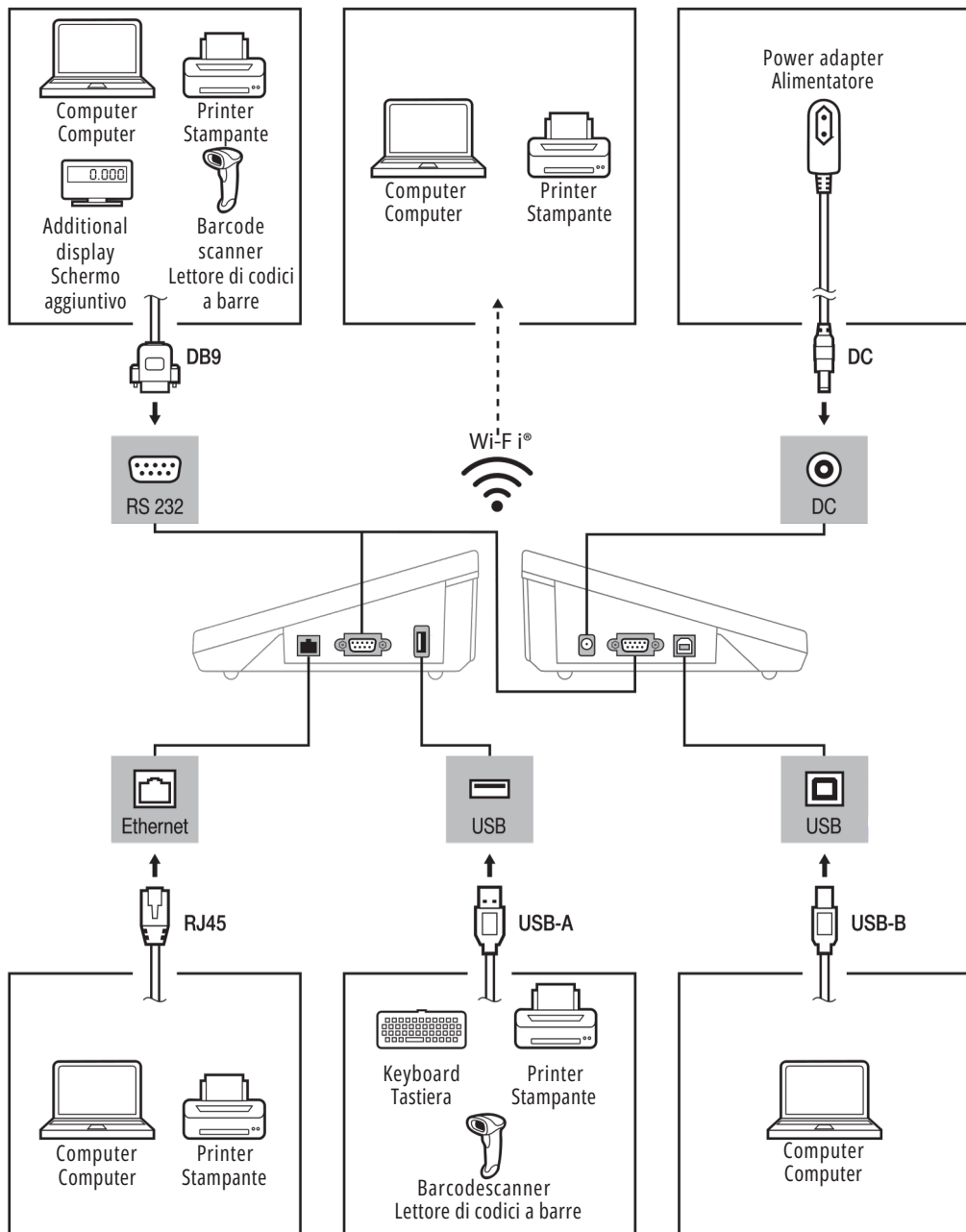
Following the above procedure you can configure sensors for other working modes.

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10. PERIPHERAL DEVICES CONNECTORS

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Wi-Fi® is a registered trademark of Wi-Fi Alliance®.

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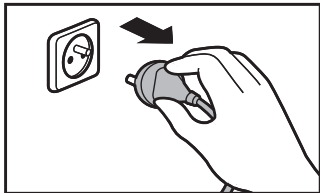
PULIZIA DEL DISPOSITIVO


11.1. General recommendations

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
Prior to cleaning it is necessary to disconnect the device from the mains! Remove the plug from the outlet, disconnect all cables.

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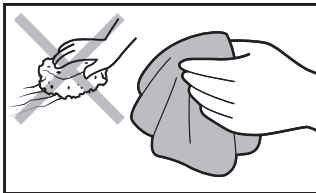Use soft cloths made of microfiber, natural fiber or man-made fiber. Avoid using abrasive cloths or cloths that might scratch the surface.

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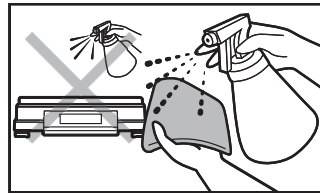Apply the cleanser onto the cloth first. Avoid applying the cleanser directly onto the device.

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Do not use aggressive cleaning agents (e.g. solvents, chlorine preparations, corrosive chemicals, bleach).

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Neither use substances of sharp structure nor intended for scrubbing.

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Make sure that neither dust nor liquid gets inside the weighing system (device inside).

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